


## Technical Information about c[G(2',5')p-2'-Fluo-AHC-A(3',5')p]

Fluorescent analogue of the metazoan cyclic dinucleotide second messenger c[G(2',5')pA(3',5')p]

Update: September 8, 2021 cw

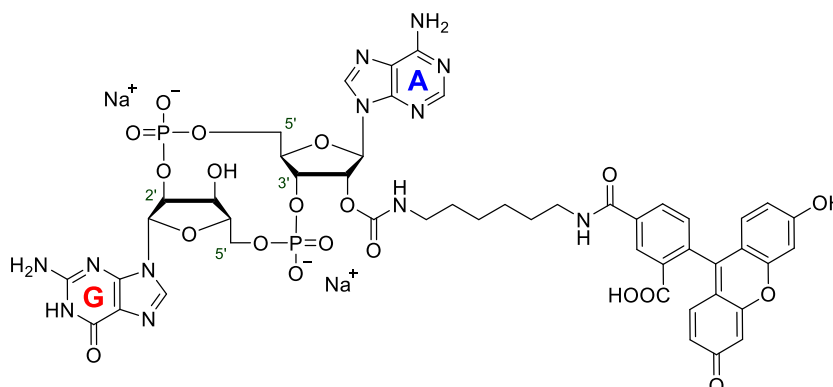

**Abbreviation:** c[G(2',5')p-2'-Fluo-AHC-A(3',5')p]

Formula	CAS No.	Molecular Weight	UV	BIOLOG Cat. No.
C <sub>48</sub> H <sub>48</sub> N <sub>12</sub> O <sub>20</sub> P <sub>2</sub> (free acid)	[pending]	1174.9 (free acid)	λ <sub>max</sub> ~494 nm / ε ~79000 / pH 9	C 195

**Name:** Cyclic (guanosine- (2' → 5')- monophosphate- 2'- O- (6- [fluoresceinyl]amino)hexylcarbamoyl)adenosine- (3' → 5')- monophosphate)

Syn.: cG-2'-Fluo-AHC-AMP(2'-5') / 2'3'-cG-2'-Fluo-AHC-AMP / 2',5'-3',5'-cG-2'-Fluo-AHC-AMP

**Description:** c[G(2',5')p-2'-Fluo-AHC-A(3',5')p] is an analogue of the metazoan cyclic dinucleotide second messenger c[G(2',5')pA(3',5')p] (2'3'-cGAMP, Cat. No. C 161) in which 5-carboxyfluorescein has been attached to the 2'-hydroxy group of the adenosine via an aminoethylcarbamoyl spacer.

**Properties:** c[G(2',5')p-2'-Fluo-AHC-A(3',5')p] is an analogue of c[G(2',5')pA(3',5')p], fluorescent with λ<sub>exc</sub> 494 nm and λ<sub>em</sub> 517 nm. The parent compound c[G(2',5')pA(3',5')p] (Cat. No. C 161) was found to be the metazoan second messenger produced by the mammalian innate immune DNA sensor cGAMP synthase (cGAS).

**Specification:** Crystallized or lyophilized sodium salt. Please keep in mind that equal concentrations of the compound may look different in volume due to sensitivity of the lyophilized form to humidity. The compound can even contract to small droplets. Normally the product is located in the conical bottom of the tube. Micromolar quantities are determined by UV/VIS at λ<sub>max</sub>.

**Purity:** Typical analysis is better than 95% (HPLC / UV 258 nm / VIS 494 nm). The product is not sterile and has not been tested for endotoxins.

**Solubility:** c[G(2',5')p-2'-Fluo-AHC-A(3',5')p] is soluble in water (≥ 1.5 mM, limits have not been determined). Please rinse tube walls carefully and preferably use ultrasonic or vortex to achieve total and uniform mixing. When opening the tube please make sure that no substance is lost within the cap.

**Stability and Storage:** c[G(2',5')p-2'-Fluo-AHC-A(3',5')p] is chemically relatively stable. Nevertheless, the compound should be protected from light and stored in the freezer (- 20° Celsius necessary, - 80° recommended), for longer storage periods preferably in freeze-dried form.

**Toxicity and Safety:** Please keep in mind, that the *in vivo* properties of this compound are not sufficiently characterized up to now. Avoid skin contact or ingestion and allow only trained personnel to handle the product. Our products are designed, developed and sold for research purposes only! They are intended for *in vitro* and nonhuman *in vivo* laboratory applications. Any other use requires approval of health authorities.

**Not for drug, household or related uses!**

### Selected References for c[G(2',5')p-2'-Fluo-AHC-A(3',5')p]:

Karanja, C.W.; Yeboah, K.S.; Ong, W.W.S.; Sintim, H.O., *RSC Chem. Biol.*, **2**, 206 - 214 (2021): "A STING-based Fluorescent Polarization Assay for Monitoring activities of Cyclic Dinucleotide Metabolizing Enzymes"

Vyskocil, S. et.al., *J.Med.Chem.*, **64**, 6902 - 6923 (2021): "Identification of Novel Carbocyclic Pyrimidine Cyclic Dinucleotide STING Agonists for Antitumor Immunotherapy Using Systemic Intravenous Route"

Miyakawa, S.; Okui, T.; Shiraishi, T.; Yoshihara, T.; Hirayama, M.; Satomi, Y.; Hamada, T.; Nishida, M.; Akimoto, C.; Sato, S.; *J. Immunol Methods*, **480**:112751 (2020): "Development of Novel Highly Sensitive Methods to Detect Endogenous cGAMP in Cells and Tissue"

Park, C.G.; Hartl, C.A.; Schmid, D.; Carmona, E.M.; Kim, H.J.; Goldberg, M.S., *Sci. Transl. Med.*, **10**(433) (2018): "Extended Release of Perioperative Immunotherapy Prevents Tumor Recurrence and Eliminates Metastases"

### Selected References for the Parent Compound c[G(2',5')pA(3',5')p]:

Lolicato, M.; Bucchi, A.; Arrigoni, C.; Zucca, S.; Nardini, M.; Schroeder, I.; Simmons, K.; Aquila, M.; DiFrancesco, D.; Bolognesi, M.; Schwede, F.; Kashin, D.; Fishwick, C.W.; Johnson, A.P.; Thiel, G.; Moroni, A., *Nat. Chem. Biol.*, **10**, 457 - 462 (2014): "Cyclic Dinucleotides bind the C-linker of HCN4 to Control Channel cAMP Responsiveness"

Konno, H.; Konno, K.; Barber, G.N., *Cell*, **155**, 688 - 698 (2013): "Cyclic Dinucleotides Trigger ULK1 (ATG1) Phosphorylation of STING to Prevent Sustained Innate Immune Signaling"

Ablasser, A.; Schmid-Burgk, J.L.; Hemmerling, I.; Horvath, G.L.; Schmidt, T.; Latz, E.; Hornung, V., *Nature*, **503**, 530 - 534 (2013): "Cell Intrinsic Immunity Spreads to Bystander Cells via the Intercellular Transfer of cGAMP"

Li, X.D.; Wu, J.; Gao, D.; Wang, H.; Sun, L.; Chen, Z.J., *Science*, **341**, 1390 - 1394 (2013): "Pivotal Roles of cGAS-cGAMP Signaling in Antiviral Defense and Immune Adjuvant Effects"

Gao, D.; Wu, J.; Wu, Y.T.; Du, F.; Aroh, C.; Yan, N.; Sun, L.; Chen, Z.J., *Science*, **341**, 903 - 906 (2013): "Cyclic GMP-AMP Synthase is an Innate Immune Sensor of HIV and Other Retroviruses"

Gao, P.; Ascano, M.; Wu, Y.; Barchet, W.; Gaffney, B.L.; Zillinger, T.; Serganov, A.A.; Liu, Y.; Jones, R.A.; Hartmann, G.; Tuschl, T.; Patel, D.J., *Cell*, **153**, 1094 - 1107 (2013): "Cyclic [G(2',5')pA(3',5')p] Is the Metazoan Second Messenger Produced by DNA-Activated Cyclic GMP-AMP Synthase"

Ablasser, A.; Goldeck, M.; Cavlar, T.; Deimling, T.; Witte, G.; Röhl, I.; Hopfner, K.-P.; Ludwig, J.; Hornung, V., *Nature*, **498**, 380 - 384 (2013): "cGAS Produces a 2'-5'-linked Cyclic Dinucleotide Second Messenger that Activates STING"

Zhang, X.; Shi, H.; Wu, J.; Zhang, X.; Sun, L.; Chen, C.; Chen, Z.J., *Mol. Cell*, **51**, 226 - 235 (2013): "Cyclic GMP-AMP Containing Mixed Phosphodiester Linkages Is An Endogenous High-Affinity Ligand for STING"

Diner, E.J.; Burdette, D.L.; Wilson, S.C.; Monroe, K.M.; Kellenberger, C.A.; Hyodo, M.; Hayakawa, Y.; Hammond, M.C.; Vance, R.E., *Cell Rep.*, **3**, 1355 - 1361 (2013): "The Innate Immune DNA Sensor cGAS Produces a Noncanonical Cyclic Dinucleotide that Activates Human STING"

Wu, J.; Sun, L.; Chen, X.; Du, F.; Shi, H.; Chen, C.; Chen, Z.J., *Science*, **339**, 826 - 830 (2013): "Cyclic GMP-AMP is an Endogenous Second Messenger in Innate Immune Signaling by Cytosolic DNA"

Sun, L.; Wu, J.; Du, F.; Chen, X.; Chen, Z.J., *Science*, **339**, 786 - 791 (2013): "Cyclic GMP-AMP Synthase is a Cytosolic DNA Sensor That Activates the Type I Interferon Pathway"