


Technical Information about Sp-2'-F-dGTP- α -S

Update: February 23, 2023ss

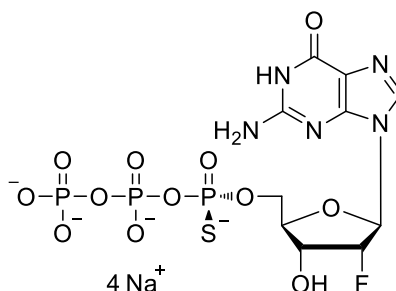

Abbreviation: **Sp-2'-F-dGTP- α -S**

Formula	CAS No.	Molecular Weight	UV	BIOLOG Cat. No.
C ₁₀ H ₁₅ FN ₅ O ₁₂ P ₃ S (free acid)	[pending]	541.2 (free acid)	λ_{\max} 252 nm / ϵ 14300 / pH 7	D 232

Name: 2'- Deoxy- 2'- fluoro guanosine - 5'- O- (1- thiotriphosphate), Sp- isomer (Sp-2'-F-dGTP- α -S), sodium salt

Description: Sp-2'-F-dGTP- α -S is an analogue of GTP and Sp-GTP- α -S, respectively. The 2'-hydroxyl group of GTP is replaced by a fluorine atom and one of the non-bridging oxygens in the S position of the α -phosphate is replaced by sulfur. The suffix "p" indicates that R/S nomenclature refers to phosphorus. The corresponding Rp-isomer is offered as well (Cat. No. D 231).

Properties:

- Increased metabolic stability compared to the parent compounds 2'-F-dGTP and GTP/dGTP, respectively,
- Useful for modulation of dGTP-responsive receptors and determination of their stereospecificity,
- May be accepted by DNA polymerases for incorporation of phosphorothioate into DNA.

Specification: Aqueous solution of the sodium salt (10 mM). Other salt forms of Sp-2'-F-dGTP- α -S are available upon request. Micromolar quantities are determined by UV at λ_{\max} . When opening the tube please make sure that no liquid is lost within the cap. A short spin-down in a bench centrifuge is recommended before use.

Purity: Typical purity is better than 95% (HPLC / UV / 252 nm) at time of quality control and packing. However, actual purity depends on storage and transport conditions. The product is not sterile and has not been tested for endotoxins.

Stability and Storage: Sp-2'-F-dGTP- α -S is most stable when stored as aqueous solution in the freezer (-20° Celsius necessary, -70° recommended), however, at ambient temperature the compound slowly starts to decompose. Thus, in order to maintain its original high quality it is recommended to allow thawing only before using the product. If you will not use up the vial with one application, please aliquot the contents of the vial in order to avoid repeated freeze/thaw cycles for the rest. When making such aliquots be sure to operate quickly and to freeze the vial again as soon as possible.

Toxicity and Safety: Since nucleoside triphosphates have important tasks in every organism, it is very likely that triphosphate analogues will interfere with many cell regulation processes *in vivo*. However, due to the rather small quantities to work with, no health hazards have been reported. Nevertheless please keep in mind, that the *in vivo* properties of this compound are not sufficiently characterized up to now. Avoid skin contact or ingestion and allow only trained personnel to handle the product.

Our products are designed, developed and sold for research purposes only! They are intended for *in vitro* and nonhuman *in vivo* laboratory applications. Any other use requires approval of health authorities.

Not for drug, household or related uses!

Reference for Sp-2'-F-dGTP- α -S: Sp-2'-F-dGTP- α -S is a new product and there are currently no corresponding references available.

Selected References for the Related Compound Sp-dGTP- α -S (Cat. No. D 031):

- Ruff, A.J.; Marienhagen, J.; Verma, R.; Roccatano, D.; Genieser, H.-G.; Nieman, P.; Shivange, A.V.; Schwaneberg, U., *J. of Mol. Cat. B: Enzymatic*, **84**, 40 - 47 (2012): "dRTP and dPTP a Complementary Nucleotide Couple for the Sequence Saturation Mutagenesis (SeSaM) Method"
- Lutz, S.; Ostermeier, M.; Benkovic, S.J., *Nucl. Acids Res.*, **29**, e16 (2001): "Rapid Generation of Incremental Truncation Libraries for Protein Engineering Using Alpha-Phosphorothioate Nucleotides"
- Sauer, S.; Lechner, D.; Berlin, K.; Lehrach, H.; Escary, J.L.; Fox, N.; Gut, I.G., *Nucleic Acids Res.*, **28** (5), e13 (2000): "A Novel Procedure for Efficient Genotyping of Single Nucleotide Polymorphisms"
- Sauer, S.; Lechner, D.; Berlin, K.; Plancon, C.; Heuermann, A.; Lehrach, H.; Gut, I.G., *Nucleic Acids Res.*, **28** (23), 100e (2000): "Full Flexibility Genotyping of Single Nucleotide Polymorphisms by the GOOD Assay"
- Xu, S.H.; Gaskin, F., *Biochim. Biophys. Acta*, **1383**, 111 - 122 (1998): "Probing the ATP Binding Site of Tubulin with Thiotriphosphate Analogues of ATP"
- Schachter, J.B.; Harden, T.K., *Br. J. Pharmacol.*, **121**, 338 - 344 (1997): "An Examination of Deoxyadenosine 5' (alpha-thio)triphosphate as a Ligand to Define P2Y Receptors and its Selectivity as a low Potency Partial Agonist of the P2Y1 Receptor"
- Nyren, P.; Karamohamed, S.; Ronaghi, M., *Anal. Biochem.*, **244**, 367 - 373 (1997): "Detection of Single-Base Changes Using a Bioluminometric Primer Extension Assay"
- Conrad, F.; Hanne, A.; Gaur, R.K.; Krupp, G., *Nucl. Acids Res.*, **23**, 1845 - 1853 (1995): "Enzymatic Synthesis of 2'-modified Nucleic Acids: Identification of Important Phosphate and Ribose Moieties in RNase P Substrates"
- Ciafre, S.A.; Rinaldi, M.; Gasparini, P.; Seripa, D.; Bisceglia, L.; Zelante, L.; Farace, M.G.; Fazio, V.M., *Nucl. Acids Res.*, **23**, 4134 - 4142 (1995): "Stability and Functional Effectiveness of Phosphorothioate Modified Duplex DNA and Synthetic "Minigenes""
- Olsen, D.B.; Sayers, J.R.; Eckstein, F., *Methods Enzymol.*, **217**, 189 - 217 (1993): "Site-directed Mutagenesis of Single-stranded and Double-stranded DNA by phosphorothioate Approach"
- King, P.; Goodburn, S., *Nucl. Acids Res.*, **20**, 1039 - 1044 (1992): "A Method for Sequence-specific Deletion Mutagenesis"
- Lee, L.G.; Connell, C.R.; Woo, S.L.; Cheng, R.D.; McArdle, B.F.; Fuller, C.W.; Halloran, N.D.; Wilson, R.K., *Nucleic Acids Res.*, **20**, 2471 - 2483 (1992): "DNA Sequencing with Dye-labeled Terminators and T7 DNA Polymerase: Effect of Dyes and dNTPs on Incorporation of Dye-terminators and Probability Analysis of Termination Fragments"
- Olsen, D.B.; Kotzorek, G.; Eckstein, F., *Biochemistry* **29**, 9546 - 9551 (1990): "Investigation of the Inhibitory Role of Phosphorothioate Internucleotidic Linkages on the Catalytic Activity of the Restriction Endonuclease EcoRV"
- Gish, G.; Eckstein, F., *Science* **240**, 1520 - 1522 (1988): "DNA and RNA Sequence Determination Based on Phosphorothioate Chemistry"
- Nakamaye, K.L.; Gish, G.; Eckstein, F.; Vosberg, H.-P., *Nucl. Acids Res.*, **16**, 9947 - 9959 (1988): "Direct Sequencing of Polymerase Chain Reaction Amplified DNA Fragments Through the Incorporation of Deoxynucleoside alpha-Thiotriphosphates"
- Abbotts, J.; SenGupta, D.N.; Zon, G.; Wilson, S.H., *J. Biol. Chem.*, **263**, 15094 - 15103 (1988): "Studies on the Mechanism of Escherichia coli DNA Polymerase I Large Fragment"
- Labeit, S.; Lehrach, H.; Goody, R.S., *Methods Enzymol.*, **155**, 166 - 177 (1987): "DNA Sequencing Using Alpha-Thiodeoxynucleotides"
- Vossberg, H.-P.; Eckstein, F., *J. Biol. Chem.*, **257**, 6595 - 6599 (1982): "Effect of Deoxynucleoside Phosphorothioates Incorporated in DNA on Cleavage by Restriction Enzymes"
- Kunkel, T.A.; Eckstein, F.; Mildvan, A.S.; Koplitz, R.M.; Loeb, L.A., *Proc. Natl. Acad. Sci. USA*, **78**, 6734 - 6738 (1981): "Deoxynucleoside [1-thio]triphosphates Prevent Proofreading During in vitro DNA Synthesis"

Selected References for the Related Compound 2'-F-dGTP:

- Richardson, F.C.; Kuchta, R.D.; Mazurkiewicz, A.; Richardson, K.A., *Biochem. Pharmacol.*, **59**, 1045 - 1052 (2000): "Polymerization of 2'-Fluoro- and 2'-O-Methyl-dNTPs by Human DNA Polymerase Alpha, Polymerase Gamma, and Primase"
- Zhu, B.; Hernandez, A.; Tan, M.; Wollenhaupt, J.; Tabor, S.; Richardson, C.C., *Nucleic Acids Res.*, **43**, e94 (2015): "Synthesis of 2'-Fluoro RNA by Syn5 RNA Polymerase"
- Ono, T.; Scalf, M.; Smith, L.M., *Nucleic Acids Res.*, **25**, 4581 - 4588 (1997): "2'-Fluoro Modified Nucleic Acids: Polymerase-Directed Synthesis, Properties and Stability to Analysis by Matrix-Assisted Laser Desorption/Ionization Mass Spectrometry"